

# Human IFN $\gamma$ High Sensitivity ELISA Kit

Instructions for use

|                    |                         |
|--------------------|-------------------------|
| Catalogue numbers: | 1x48 tests: 850.900.048 |
|                    | 1x96 tests: 850.900.096 |
|                    | 2x96 tests: 850.900.192 |

**For research use only**

Fast Track Your Research.....

## Table of Contents

|                                                              |    |
|--------------------------------------------------------------|----|
| 1. Intended use.....                                         | 2  |
| 2. Principle of the method.....                              | 2  |
| 3. Reagents provided and reconstitution .....                | 2  |
| 4. Materials required but not provided .....                 | 3  |
| 5. Storage Instructions.....                                 | 3  |
| 6. Specimen collection, processing & storage .....           | 3  |
| 7. Safety & precautions for use .....                        | 4  |
| 8. Assay Preparation .....                                   | 5  |
| 8.1. Assay Design.....                                       | 5  |
| 8.2. Preparation of Wash Buffer .....                        | 5  |
| 8.3. Preparation of Standard Diluent Buffer.....             | 5  |
| 8.4. Preparation of Standard .....                           | 6  |
| 8.5. Preparation of Samples .....                            | 6  |
| 8.6. Preparation of Biotinylated anti-IFN $\gamma$ .....     | 6  |
| 8.7. Preparation of Streptavidin-HRP solutions 1 and 2 ..... | 7  |
| 8.8. Preparation of Amplifier .....                          | 7  |
| 9. Method.....                                               | 8  |
| 10. Data Analysis.....                                       | 9  |
| 11. Assay limitations .....                                  | 10 |
| 12. Performance Characteristics .....                        | 10 |
| 12.1. Sensitivity.....                                       | 10 |
| 12.2. Specificity .....                                      | 10 |
| 12.3. Precision.....                                         | 10 |
| 12.4. Dilution Parallelism .....                             | 10 |
| 12.5. Spike Recovery.....                                    | 11 |
| 12.6. Expected Values.....                                   | 11 |
| 13. Assay Summary.....                                       | 12 |

# Human IFN $\gamma$ High Sensitivity ELISA KIT

## 1. Intended use

The Human high sensitivity IFN $\gamma$  ELISA Kit is to be used for the in-vitro quantitative determination of interferon gamma (IFN $\gamma$ ) in human sera, plasmas, buffered solutions or cell culture media. The assay will recognize both natural and recombinant human IFN $\gamma$ .

**This kit has been configured for research use only.**

## 2. Principle of the method

The IFN $\gamma$  Kit is a solid phase sandwich Enzyme Linked-Immuno- Sorbent Assay (ELISA). A monoclonal antibody specific for IFN $\gamma$  has been coated onto the wells of the microtiter strips provided. Samples, including standards of known IFN $\gamma$  concentrations and unknowns are pipetted into these wells.

During the first incubation, the IFN $\gamma$  antigen and a biotinylated monoclonal antibody specific for IFN $\gamma$  are successively incubated.

After washing, the enzyme (streptavidin-peroxydase) is added. All the unbound enzyme is removed by washing and the first amplification step is performed by adding the Biotine-Tyramine reagent. Under the action of HRP, a biotine polymerisation reaction occurs in the region of the HRP linked to the detection antibody. After washing the second amplification step is performed and the polymerised biotine is revealed by a new streptavidin-HRP step. Finally after washing, the substrate is added. The intensity of this coloured product is directly proportional to the concentration of IFN $\gamma$  present in the samples.

## 3. Reagents provided and reconstitution

| Reagents<br>(Store@2-8°C)                   | Quantity<br>1x48 well kit<br>Cat no.<br>850.900.048 | Quantity<br>1x96 well kit<br>Cat no.<br>850.900.096 | Quantity<br>2x96 well kit<br>Cat no.<br>850.900.192 | Reconstitution                                                                     |
|---------------------------------------------|-----------------------------------------------------|-----------------------------------------------------|-----------------------------------------------------|------------------------------------------------------------------------------------|
| 96 well microtitre strip plate              | 1/2                                                 | 1                                                   | 2                                                   | Ready to use (Pre-coated)                                                          |
| Plastic plate covers                        | 2                                                   | 2                                                   | 4                                                   | n/a                                                                                |
| Standard: 25 pg/ml                          | 1                                                   | 2                                                   | 4                                                   | Reconstitute as directed on the vial<br>(see Assay preparation, section 8)         |
| Standard Diluent (Buffer)                   | 1 (25ml)                                            | 1 (25ml)                                            | 1 (25ml)                                            | 10x Concentrate, dilute in distilled water<br>(see reagent preparation, section 8) |
| Biotinylated anti-IFN $\gamma$              | 1 (0.4ml)                                           | 1 (0.4ml)                                           | 2 (0.4ml)                                           | Dilute in Biotinylated Antibody Diluent<br>(see Assay preparation, section 8)      |
| Biotinylated Antibody Diluent               | 1 (7.5ml)                                           | 1 (7.5ml)                                           | 1 (13ml)                                            | Ready to use                                                                       |
| Streptavidin-HRP                            | 1 (5 $\mu$ l)                                       | 2 (5 $\mu$ l)                                       | 4 (5 $\mu$ l)                                       | Add 0.5ml of HRP diluent prior to use<br>(see Assay preparation, section 8)        |
| Amplification Diluent                       | 1 (25ml)                                            | 1 (25ml)                                            | 1 (25ml)                                            | Ready to use                                                                       |
| Amplifier*                                  | 1 (200 $\mu$ l)                                     | 1 (200 $\mu$ l)                                     | 2 (200 $\mu$ l)                                     | Dilute in Amplification Buffer                                                     |
| HRP Diluent                                 | 1 (25ml)                                            | 1 (25ml)                                            | 2 (25ml)                                            | Ready to use                                                                       |
| Wash Buffer                                 | 1 (10ml)                                            | 1 (10ml)                                            | 2 (10ml)                                            | 200x Concentrate dilute in distilled water<br>(see Assay preparation, section 8)   |
| TMB Substrate                               | 1 (11ml)                                            | 1 (11ml)                                            | 1 (24ml)                                            | Ready to use                                                                       |
| H <sub>2</sub> SO <sub>4</sub> stop reagent | 1 (11ml)                                            | 1 (11ml)                                            | 2 (11ml)                                            | Ready to use                                                                       |

\*Reagent contains ethyl alcohol

## 4. Materials required but not provided

- Microtiter plate reader fitted with appropriate filters (450nm required with optional 630nm reference filter)
- Microplate washer or wash bottle
- 10, 50, 100, 200 and 1,000µl adjustable single channel micropipettes with disposable tips
- 50-300µl multi-channel micropipette with disposable tips
- Multichannel micropipette reagent reservoirs
- Distilled water
- Vortex mixer
- Orbital shaker
- Miscellaneous laboratory plastic and/or glass, if possible sterile

## 5. Storage Instructions

Store kit reagents between 2 and 8°C. Immediately after use remaining reagents should be returned to cold storage (2-8°C). Expiry of the kit and reagents is stated on box front labels. The expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, the reagent is not contaminated by the first handling.

**Wash Buffer:** Once prepared store at 2-8°C for up to 1 week

**Standard Diluent Buffer:** Once prepared store at 2-8°C for up to 1 week

**Standards :** Once prepared use immediately and do not store

**Biotinylated Secondary Antibody:** Once prepared use immediately and do not store

**Streptavidin-HRP:** Once prepared use immediately and do not store

## 6. Specimen collection, processing & storage

Cell culture supernatants, serum, plasma or other biological samples will be suitable for use in the assay. Remove serum from the clot or red cells, respectively, as soon as possible after clotting and separation.

**Cell culture supernatants:** Remove particulates and aggregates by spinning at approximately 1000 x g for 10 min.

**Serum:** Avoid any unintentional stimulation of the cells by the procedure. Use pyrogen/endotoxin free collecting tubes. Serum should be removed rapidly and carefully from the red cells after clotting. For that, after clotting, centrifuge at approximately 1000 x g for 10 min and remove serum.

**Plasma:** EDTA, citrate and heparin plasma can be assayed. Spin samples at 1000 x g for 30 min to remove particulates. Harvest plasma.

**Storage:** If not analyzed shortly after collection, samples should be aliquoted (250-500µl) to avoid repeated freeze-thaw cycles and stored frozen at -70°C. Avoid multiple freeze-thaw cycles of frozen specimens.

**Recommendation:** Do not thaw by heating at 37°C or 56°C. Thaw at room temperature and make sure that sample is completely thawed and homogeneous before use. When possible avoid use of badly haemolysed or lipemic sera. If large amounts of particles are present these should be removed prior to use by centrifugation or filtration.

## 7. Safety & precautions for use

- Handling of reagents, serum or plasma specimens should be in accordance with local safety procedures , e.g.CDC/NIH Health manual : " Biosafety in Microbiological and Biomedical Laboratories" 1984
- Laboratory gloves should be worn at all times
- Avoid any skin contact with H<sub>2</sub>SO<sub>4</sub> and TMB. In case of contact, wash thoroughly with water
- Do not eat, drink, smoke or apply cosmetics where kit reagents are used
- Do not pipette by mouth
- When not in use, kit components should be stored refrigerated or frozen as indicated on vials or bottles labels
- All reagents should be warmed to room temperature before use. Lyophilized standards should be discarded after use
- Once the desired number of strips has been removed, immediately reseal the bag to protect the remaining strips from deterioration
- Cover or cap all reagents when not in use
- Do not mix or interchange reagents between different lots
- Do not use reagents beyond the expiration date of the kit
- Use a clean disposable plastic pipette tip for each reagent, standard, or specimen addition in order to avoid cross contamination, for the dispensing of H<sub>2</sub>SO<sub>4</sub> and substrate solution, avoid pipettes with metal parts
- Use a clean plastic container to prepare the washing solution
- Thoroughly mix the reagents and samples before use by agitation or swirling
- All residual washing liquid must be drained from the wells by efficient aspiration or by decantation followed by tapping the plate forcefully on absorbent paper. Never insert absorbent paper directly into the wells
- The TMB solution is light sensitive. Avoid prolonged exposure to light. Also, avoid contact of the TMB solution with metal to prevent colour development. Warning TMB is toxic avoid direct contact with hands. Dispose off properly
- If a dark blue colour develops within a few minutes after preparation, this indicates that the TMB solution has been contaminated and must be discarded. Read absorbance's within 1 hour after completion of the assay
- When pipetting reagents, maintain a consistent order of addition from well-to-well. This will ensure equal incubation times for all wells
- Follow incubation times described in the assay procedure
- Dispense the TMB solution within 15 min of the washing of the microtitre plate

## 8. Assay Preparation

Bring all reagents to room temperature before use

### 8.1. Assay Design

Determine the number of microwell strips required to test the desired number of samples plus appropriate number of wells needed for running zeros and standards. Each sample, standard and zero should be tested **in duplicate**. Remove sufficient Microwell Strips for testing from the pouch immediately prior to use. Return any wells not required for this assay with desiccant to the pouch. Seal tightly and return to 2-8°C storage.

**Example plate layout** (example shown for a 6 point standard curve)

|   | Standards<br>(pg/mL) |      | Sample Wells |   |   |   |   |   |   |    |    |    |
|---|----------------------|------|--------------|---|---|---|---|---|---|----|----|----|
|   | 1                    | 2    | 3            | 4 | 5 | 6 | 7 | 8 | 9 | 10 | 11 | 12 |
| A | 25                   | 25   |              |   |   |   |   |   |   |    |    |    |
| B | 12.5                 | 12.5 |              |   |   |   |   |   |   |    |    |    |
| C | 6.25                 | 6.25 |              |   |   |   |   |   |   |    |    |    |
| D | 3.12                 | 3.12 |              |   |   |   |   |   |   |    |    |    |
| E | 1.56                 | 1.56 |              |   |   |   |   |   |   |    |    |    |
| F | 0.78                 | 0.78 |              |   |   |   |   |   |   |    |    |    |
| G | zero                 | zero |              |   |   |   |   |   |   |    |    |    |
| H |                      |      |              |   |   |   |   |   |   |    |    |    |

*All remaining empty wells can be used to test samples in duplicate*

### 8.2. Preparation of Wash Buffer

Dilute the (200x) wash buffer concentrate 200 fold with distilled water to give a 1x working solution. Pour entire contents (10 ml) of the Washing Buffer Concentrate into a clean 2,000 ml graduated cylinder. Bring final volume to 2,000 ml with glass-distilled or deionized water. Mix gently to avoid foaming. Transfer to a clean wash bottle and store at 2°-8°C for up to 1 week.

### 8.3. Preparation of Standard Diluent Buffer

Add the contents of the vial (10x concentrate) to 225ml of distilled water before use.

This solution can be stored at 2-8°C for up to 1 week.

## 8.4. Preparation of Standard

Standard vials must be reconstituted with the volume of standard diluent shown on the vial immediately prior to use. This reconstitution gives a stock solution of 25 pg/ml of IFN $\gamma$ . **Mix the reconstituted standard gently by inversion only.** Serial dilutions of the standard are made directly in the assay plate to provide the concentration range from 25 to 0.78 pg/ml. A fresh standard curve should be produced for each new assay.

- Immediately after reconstitution add 200 $\mu$ l of the reconstituted standard to wells A1 and A2, which provides the highest concentration standard at 25 pg/ml
- Add 100 $\mu$ l of Standard Diluent to the remaining standard wells B1 and B2 to F1 and F2
- Transfer 100 $\mu$ l from wells A1 and A2 to B1 and B2. Mix the well contents by repeated aspirations and ejections taking care not to scratch the inner surface of the wells
- Continue this 1:1 dilution using 100 $\mu$ l from wells B1 and B2 through to wells F1 and F2 providing a serial diluted standard curve ranging from 25 pg/ml to 0.78 pg/ml
- Discard 100 $\mu$ l from the final wells of the standard curve (F1 and F2)

Alternatively these dilutions can be performed in separate clean tubes and immediately transferred directly into the relevant wells.

## 8.5. Preparation of Samples

Normal sera and plasmas may be applied undiluted. Nevertheless, sera or plasmas from patients with various pathologies may be applied undiluted and diluted (to prevent too high concentrations). Please note that certain sera/plasma may induce false positive (for example, in reason to presence of anti-mouse anti-IgG antibody). A simple sample dilution (1:2) allows eliminating interference. As IFN $\gamma$  concentrations may vary considerably in cell supernatant samples, it is not easy to recommend a dilution factor.

## 8.6. Preparation of Biotinylated anti-IFN $\gamma$

It is recommended this reagent is prepared immediately before use. Dilute the biotinylated anti- IFN $\gamma$  with the biotinylated antibody diluent in an appropriate clean glass vial using volumes appropriate to the number of required wells. Please see example volumes below:

| Number of wells required | Biotinylated Antibody ( $\mu$ l) | Biotinylated Antibody Diluent ( $\mu$ l) |
|--------------------------|----------------------------------|------------------------------------------|
| 16                       | 40                               | 1060                                     |
| 24                       | 60                               | 1590                                     |
| 32                       | 80                               | 2120                                     |
| 48                       | 120                              | 3180                                     |
| 96                       | 240                              | 6360                                     |

## 8.7. Preparation of Streptavidin-HRP solutions 1 and 2

It is recommended to centrifuge vial for a few seconds in a microcentrifuge to collect all the volume at the bottom.

Dilute the 5 $\mu$ l vial with 0.5ml of HRP diluent **immediately before use**. This pre-dilution will be used for Step 8 and step 13. Do-not keep this diluted vial for future experiments. Further dilute the HRP solution to volumes appropriate for the number of required wells in a clean glass vial. Please see example volumes below:

| Number of wells required | Streptavidin-HRP solution 1 (Step 8) |                               | Streptavidin-HRP solution 2 (Step 13) |                               |
|--------------------------|--------------------------------------|-------------------------------|---------------------------------------|-------------------------------|
|                          | Streptavidin-HRP ( $\mu$ l)          | Streptavidin-HRP Diluent (ml) | Streptavidin-HRP ( $\mu$ l)           | Streptavidin-HRP Diluent (ml) |
| 16                       | 20                                   | 1.980                         | 40                                    | 1.960                         |
| 32                       | 40                                   | 3.960                         | 80                                    | 3.920                         |
| 48                       | 60                                   | 5.940                         | 120                                   | 5.880                         |
| 96                       | 120                                  | 11.880                        | 240                                   | 11.760                        |

## 8.8. Preparation of Amplifier

It is recommended this reagent is prepared immediately before use. Dilute the Amplifier with the Amplification diluent in an appropriate clean glass vial using volumes appropriate to the number of required wells. Please see example volumes below:

| Number of wells required | Amplifier ( $\mu$ l) | Amplification Diluent (ml) |
|--------------------------|----------------------|----------------------------|
| 16                       | 20                   | 1.980                      |
| 32                       | 40                   | 3.960                      |
| 48                       | 60                   | 5.940                      |
| 96                       | 120                  | 11.880                     |

## 9. Method

We strongly recommend that every vial is mixed thoroughly without foaming prior to use except the standard vial which must be mixed gently by inversion only. Prepare all reagents as shown in section 8.

**Note:** Final preparation of Biotinylated anti-IFN $\gamma$  (section 8.6) and Streptavidin-HRP (section 8.7) should occur immediately before use.

| Assay Step                                                                                                                                                                                                                    |            | Details                                                                                                                                                                                                                                          |
|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| 1.                                                                                                                                                                                                                            | Addition   | Add 100 $\mu$ l of each <b>standard, sample and zero</b> (Standard Diluent Buffer) in duplicate to appropriate number of wells                                                                                                                   |
| 2.                                                                                                                                                                                                                            | Incubation | Cover with a plastic plate cover and incubate at room temperature (18 to 25°C) with slow shaking for <b>1 hour</b>                                                                                                                               |
| 3.                                                                                                                                                                                                                            | Wash       | Remove the cover and wash the plate as follows:<br>a) Aspirate the liquid from each well<br>b) Dispense 0.3 ml of <b>1x washing solution</b> into each well<br>c) Aspirate the contents of each well<br>d) Repeat step b and c another two times |
| 4.                                                                                                                                                                                                                            | Addition   | Add 50 $\mu$ l of diluted <b>biotinylated anti-IFN<math>\gamma</math></b> to all wells                                                                                                                                                           |
| 5.                                                                                                                                                                                                                            | Incubation | Cover with a plastic plate cover and incubate at room temperature (18 to 25°C) with slow shaking for <b>1 hour</b>                                                                                                                               |
| 6.                                                                                                                                                                                                                            | Wash       | Repeat wash step 3.                                                                                                                                                                                                                              |
| 7.                                                                                                                                                                                                                            | Addition   | Add 100 $\mu$ l of <b>Streptavidin-HRP solution 1</b> into all wells                                                                                                                                                                             |
| 8.                                                                                                                                                                                                                            | Incubation | Cover with a plastic plate cover and incubate at room temperature (18 to 25°C) with slow shaking for <b>20 min</b>                                                                                                                               |
| 9.                                                                                                                                                                                                                            | Wash       | Repeat wash step 3.                                                                                                                                                                                                                              |
| 10.                                                                                                                                                                                                                           | Addition   | Add 100 $\mu$ l of diluted <b>Amplifier</b> to all wells                                                                                                                                                                                         |
| 11.                                                                                                                                                                                                                           | Incubation | Cover with a plastic plate cover and incubate at room temperature (18 to 25°C) with slow shaking for <b>15 min</b>                                                                                                                               |
| 12.                                                                                                                                                                                                                           | Wash       | Repeat wash step 3.                                                                                                                                                                                                                              |
| 13.                                                                                                                                                                                                                           | Addition   | Add 100 $\mu$ l of <b>Streptavidin-HRP solution 2</b> into all wells                                                                                                                                                                             |
| 14.                                                                                                                                                                                                                           | Incubation | Cover with a plastic plate cover and incubate at room temperature (18 to 25°C) with slow shaking for <b>20 min</b>                                                                                                                               |
| 15.                                                                                                                                                                                                                           | Wash       | Repeat wash step 3.                                                                                                                                                                                                                              |
| 16.                                                                                                                                                                                                                           | Addition   | Add 100 $\mu$ l of ready-to-use <b>TMB Substrate Solution</b> into all wells                                                                                                                                                                     |
| 17.                                                                                                                                                                                                                           | Incubation | Incubate in the dark for <b>10-20 minutes*</b> at room temperature. Avoid direct exposure to light by wrapping the plate in aluminium foil                                                                                                       |
| 18.                                                                                                                                                                                                                           | Addition   | Add 100 $\mu$ l of <b>H<sub>2</sub>SO<sub>4</sub>:Stop Reagent</b> into all wells                                                                                                                                                                |
| <b>Read the absorbance</b> value of each well (immediately after step 11.) on a spectrophotometer using 450 nm as the primary wavelength and optionally 630 nm as the reference wave length (610 nm to 650 nm is acceptable). |            |                                                                                                                                                                                                                                                  |

*\*Incubation time of the substrate solution is usually determined by the ELISA reader performance. Many ELISA readers only record absorbance up to 2.0 O.D. Therefore the colour development within individual microwells must be observed by the analyst, and the substrate reaction stopped before positive wells are no longer within recordable range*

**Note:** In case of incubation without shaking the O.D values may be lower than with shaking; in this case let the color develop longer in order to obtain correct OD values

## 10. Data Analysis


Calculate the average absorbance values for each set of duplicate standards and samples. Ideally duplicates should be within 20% of the mean.

Generate a linear standard curve by plotting the average absorbance of each standard on the vertical axis versus the corresponding Human IFN $\gamma$  standard concentration on the horizontal axis.

The amount of IFN $\gamma$  in each sample is determined by extrapolating OD values against IFN $\gamma$  standard concentrations using the standard curve.

**Example IFN $\gamma$  Standard curve**

| Standard | IFN $\gamma$ Conc<br>pg/ml | OD<br>(450nm)<br>Mean | CV (%) |
|----------|----------------------------|-----------------------|--------|
| 1        | 25                         | 2.297                 | 2.1    |
| 2        | 12.5                       | 1.172                 | 0.2    |
| 3        | 6.25                       | 0.712                 | 3.9    |
| 4        | 3.12                       | 0.487                 | 3.1    |
| 5        | 1.56                       | 0.381                 | 2.4    |
| 6        | 0.78                       | 0.322                 | 5.3    |
| Zero     | 0                          | 0.236                 | 7.8    |

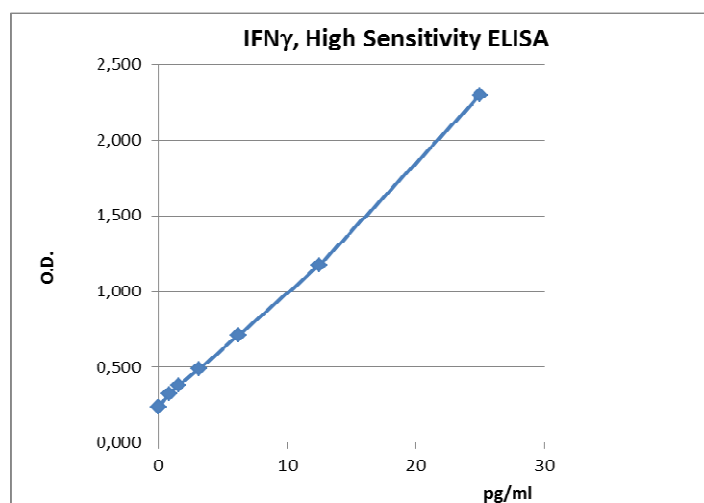

**Note:** curve shown above should not be used to determine results. Every laboratory must produce a standard curve for each set of microwell strips assayed.

## 11. Assay limitations

Do not extrapolate the standard curve beyond the maximum standard curve point. The dose-response is non-linear in this region and good accuracy is difficult to obtain. Concentrated samples above the maximum standard concentration must be diluted with Standard diluent or with your own sample buffer to produce an OD value within the range of the standard curve. Following analysis of such samples always multiply results by the appropriate dilution factor to produce actual final concentration.

The influence of various drugs on end results has not been investigated. Bacterial or fungal contamination and laboratory cross-contamination may also cause irregular results.

Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Completely empty wells before dispensing fresh Washing Buffer, fill with Washing Buffer as indicated for each wash cycle and do not allow wells to sit uncovered or dry for extended periods.

Disposable pipette tips, flasks or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.

As with most biological assays conditions may vary from assay to assay therefore **a fresh standard curve must be prepared and run for every assay.**

## 12. Performance Characteristics

### 12.1. Sensitivity

The sensitivity, minimum detectable dose of Human IFN $\gamma$  using this Diaclone Human IFN $\gamma$  High Sensitivity ELISA kit was found to be **<0.69 pg/ml**. This was determined by adding 3 standard deviations to the mean OD obtained when the zero standard was assayed 40 times.

### 12.2. Specificity

Ten specificities were tested with concentrations higher than IFN $\gamma$  curve concentrations. No cross reaction was observed for concentrations ranging from 2500 to 78.12 pg/ml for IL-1 $\alpha$ , IL-2, IL-8, IL-12p40, TNF $\alpha$ , CD95 Fas, TRAIL, ICAM-1, Gp130 and GM-CSF.

### 12.3. Precision

Two human serum pools, two human plasma pools and two cell culture media samples with various concentrations of IFN $\gamma$  were tested for repeatability and reproductibility. Each assay was carried out with 3 duplicates of each sample. Three independent assays were performed. The intra-assay and inter-assay coefficient of variation has been calculated to be 3.9% and 8.6% respectively.

### 12.4. Dilution Parallelism

Two human serum pools, one human plasma pool and one cell culture medium samples with various concentrations of IFN $\gamma$  were serially diluted in standard buffer diluent. Linearity was evaluated on 4 dilutions. The linear regression of samples versus the expected concentrations yielded a quote slope of 0.994.

## **12.5. Spike Recovery**

The spike recovery was evaluated by spiking two levels of IFN $\gamma$  into three human serum pools, two plasma pools and two culture media. Recovery was evaluated with one test. Mean recoveries were 92% in sera (range values: 84-98%), 72% in plasma (Range values: 61-81%) and 108% in cell culture media (range values: 88-117%).

## **12.6. Expected Values**

16 sera and 16 plasmas from healthy individual donors were tested undiluted and diluted in duplicates. Mean concentration of IFN $\gamma$  in sera was 2.3 pg/ml (12 positives – range values of positive sera : 0.70 – 5.82 pg/ml) and in plasma is 1.89 pg/ml (13 positives – range values of positive plasmas : 0.89 – 5.08 pg/ml).

## 13. Assay Summary

Total procedure length : 3h00mn

|                                         |                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                        |
|-----------------------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Specific molecule detection steps       | <p>Add 100µl of sample or diluted standard</p> <p>↓</p> <p>Incubate 1 hour at room temperature with slow shaking</p> <p>↓</p> <p>Wash three times</p> <p>↓</p> <p>Add 50µl of diluted biotinylated detection antibody to all wells</p> <p>↓</p> <p>Incubate 1 hour at room temperature with slow shaking</p>                                                                                                                                                                                                           |
| Streptavidin –HRP / Amplification steps | <p>↓</p> <p>Wash three times</p> <p>↓</p> <p>Add 100µl of streptavidin-HRP (solution 1) to all wells</p> <p>↓</p> <p>Incubate 20 min at room temperature with slow shaking</p> <p>↓</p> <p>Wash three times</p> <p>↓</p> <p>Add 100 µl amplifier to all wells</p> <p>↓</p> <p>Incubate 15 min at room temperature with slow shaking</p> <p>↓</p> <p>Wash three times</p> <p>↓</p> <p>Add 100µl of streptavidin-HRP (solution 2) to all wells</p> <p>↓</p> <p>Incubate 20 min at room temperature with slow shaking</p> |
| Revelation and reading steps            | <p>↓</p> <p>Wash three times</p> <p>↓</p> <p>Add 100 µl of ready-to-use TMB</p> <p>Protect from light. Let the color develop for around 5 min.</p> <p>↓</p> <p>Add 100 H<sub>2</sub>SO<sub>4</sub></p> <p>↓</p> <p>Read Absorbance at 450 nm</p>                                                                                                                                                                                                                                                                       |

### Products Manufactured and Distributed by:

Diaclone SAS  
 1 Boulevard A.Fleming  
 25020 Besancon Cedex  
 France  
 Tel +33 381413838  
 Fax +33 381413636  
 Email: [info@diaclone.com](mailto:info@diaclone.com)